#### related products:

Cat. No. CP-0110

CentriPure P2
Gel filtration column

for 200 µl sample volume

Cat. No. CP-0107

**CentriPure P10** 

Gel filtration column for 1.0 ml sample volume

Cat. No. CP-0108

CentriPure P25
Gel filtration column

for 2.5 ml sample volume

Cat. No. CP-0113

**CentriPure P50** 

Gel filtration column for 5.0 ml sample volume

Cat. No. CP-0119

CentriPure P100

Gel filtration column for 10 ml sample volume

Cat. No. CP-0131

CentriPure P500

Gel filtration column for 50 ml sample volume

Cat. No. CP-0132

**CentriPure Dolly Mix** 

Assorted columns for protein purification
Five each of P2, P5, P10 and P25 columns



LabRack for CentriPure columns

The LabRack column processing station makes purification easy and convenient



emp BIOTECH GmbH · Robert-Rössle-Str. 10 · 13125 Berlin · Germany Tel. +49 (0)30 94 89 22 01 · info@empbiotech.com · www.empbiotech.com

emp BIOTECH GmbH is an ISO 9001:2008 certified company registration number 011001300789 (TÜV Rheinland)



Cat. No. CP-0106

#### **CentriPure P5**

Hydrated gel filtration column for protein purification



Instructions for use

## 1. Column Preparation



Remove the cap from the top and then the bottom cap of the **CentriPure P5** column.


Allow excess column fluid to drain (via gravity) into a suitable waste reservoir.

### 2. Column Equilibration



To equilibrate the column, allow the equilibration buffer to enter the gel bed completely and continue elution until approximately 10 ml of buffer has been eluted.

# 3. Sample Application

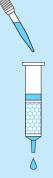

Transfer 0.5 ml sample to the **CentriPure P5** column.


Allow the sample to enter the gel bed completely.

### 4. Elution




Place a tube
for sample
collection under the
CentriPure P5 column.
Transfer 1 ml of elution
buffer to the column and
elute the purified sample.

### typical applications:

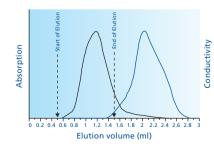

Desalting of protein solution (0.5 mg albumin (OvA) in 0.5 ml 0.8 M NaCl), elution with water

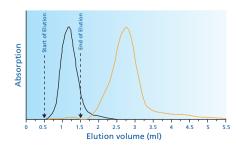

Elution profile overlay of albumin (0.5 mg OvA, black line – 280 nm) and free dye (0.25 µmol FAM, orange line – 490 nm) in NaHCO<sub>3</sub>, elution with water (0.5 ml sample volume)